pure clean origin

liquid handling

/// 2012

made in germany  made by ritter

Ripette pro and Ritips professional

Ripette genx and Ritips professional

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clinical products

About realisation and idea

For more than 20 years Ritter has been developing consumables and instruments which make liquid handling processes such as serial dispensing more convenient. In this respect our Ritter dispensers have been constructed in line with our consumables in order to continuously achieve the necessary precision.

Therefore numerous laboratories with sensible instruments are equipped with corresponding disposables from Ritter Medical, but also customer demands for new products will be developed with competence and economic efficiency, from the initial idea up to mass production.

Ritter's success is based on the independence of a medium-sized, German and family-owned company with high quality standards. All these factors make Ritter Medical a reliable partner for your requirements. Convince yourself of the new product range and the possibilities for constructive developments offered by Ritter Medical and do not hesitate to benefit of innovations developed by Ritter.

We would be pleased to provide you with competent advise.

Michael Fuchs,
Teamleader, Ritter Medical

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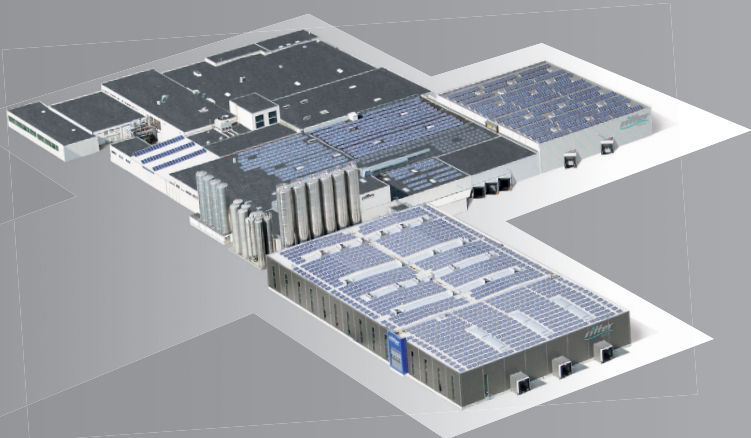
Ecological | Ritter GmbH only applies certified, pollution-free plastics. When applicable, biodegradable as well as recyclable plastics are used for the production of high-quality clinical products from Ritter.

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Product development | *Ritter GmbH was founded in 1965. The medium-sized independent family-owned company is based in Schwabmünchen, Bavaria.*

At its 25,000 m² production facility the company's 280 employees develop and produce high quality plastic products for the world market certified under DIN EN ISO 9001:2000 and 13485:2003.

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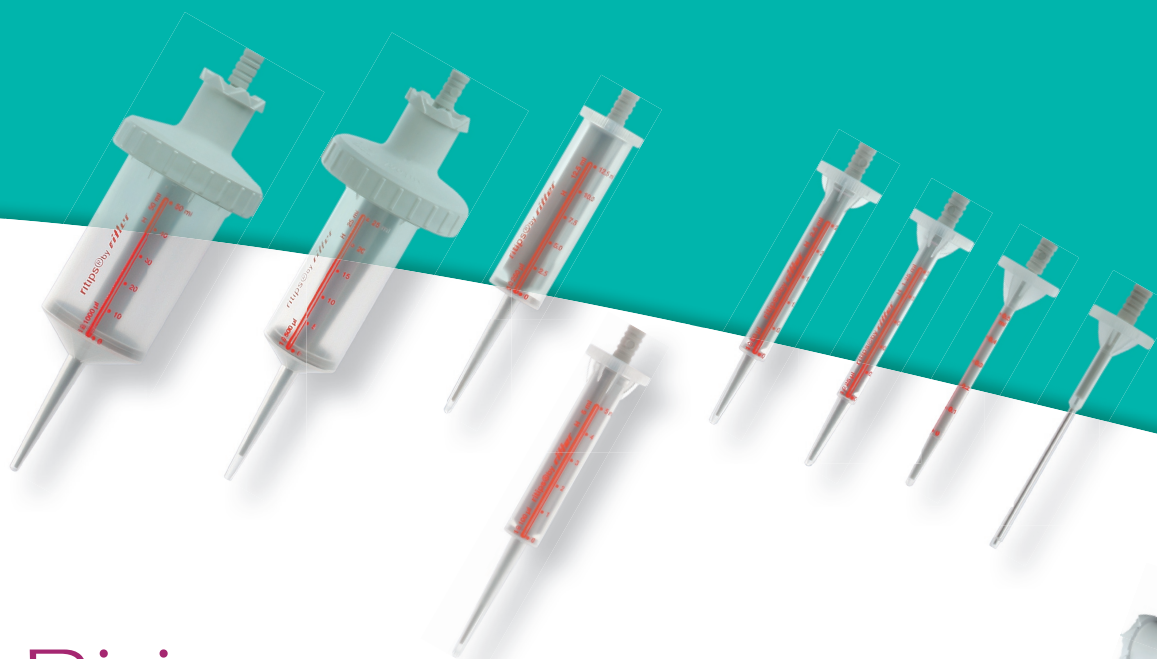
*For routine work Ritter offers the manual dispenser **Ripette**, a robust and time-saving device for long dispensing series. With this stepper up to 40 different dosing volumes can be chosen when using in total 5 different settings. With the respective **Ritips** that are available in 8 different tips sizes the optimal and cost-saving consumables can be used.*


***Ripette pro** is currently the lightest mechanical stepper worldwide with an outstanding ergonomics – both parameters are mandatory for a fatigue-free dispensing of long series. For in total 12 different available tips sizes as from 1 µl up to 5.500 µl more than 120 dosing volumes can be adjusted.*

*Thanks to the electronic support of the **Ripette genX** more than 2.500 different volumes can be dispensed with a maximum in accuracy.*

*With the **Ritips professional** Ritter offers a new generation of dispenser tips in 9 different sizes which meet the demands in research or diagnostics.*

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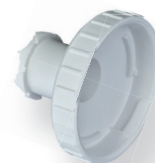Flexible: With the Ripette pro you can work flexible and practical with 12 tip sizes, 10 volume adjustments and 120 program steps.

Clear: Volume adjustments and dosing volumes for different tips are shown in a vision panel on the Ripette pro.

Handy: A perfect ergonomic design makes an easy and comfortable use possible. With the easy-moving dosing button in combination with the low weight, long dosing series can be dispensed conveniently.

Compatible: The Ripette pro is compatible not only with PD-tips from Brand but also with combitips and combitips plus from Eppendorf and has been precisely aligned to Ritips professional from Ritter.

Independent: In addition to the handy design, the Ripette pro has been constructed consistent mechanical. The Ripette pro is maintenance-free and does not need any additional resources such as power supply or batteries.

Tip size	Program number									
	1	2	3	4	5	6	7	8	9	10
0,05 ml	1	1,5	2	2,5	3	3,5	4	4,5	5	5,5
0,10 ml	2	3	4	5	6	7	8	9	10	11
0,20 ml	4	6	8	10	12	14	16	18	20	22
0,50 ml	10	15	20	25	30	35	40	45	50	55
1,00 ml	20	30	40	50	60	70	80	90	100	110
1,25 ml	25	37,5	50	62,5	75	87,5	100	112,5	125	137,5
2,50 ml	50	75	100	125	150	175	200	225	250	275
5,00 ml	100	150	200	250	300	350	400	450	500	550
10,00 ml	200	300	400	500	600	700	800	900	1000	1100
12,50 ml	250	375	500	625	750	875	1000	1125	1250	1375
25,00 ml	500	750	1000	1250	1500	1750	2000	2250	2500	2750
50,00 ml	1000	1500	2000	2500	3000	3500	4000	4500	5000	5500
	48	35	23	19	15	13	11	9	8	7
	Steps									

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Contamination-free: Ripette pro and Ritips

The Ripette pro is predestinated for dispensing of contaminated fluids and volatile liquids. Due to the exclusive acceptance of dispenser tips with plunger e.g. Ritips and other compatible tips, the Ripette pro is secure from contamination by ascending aerosols in the tip and guarantees a long-term secure use even with samples of high fluidity.

Artikel-Nr.	Artikelbezeichnung	VE	Stück / Umkarton
Article-no.	Article name	PU	Pieces/ cover. box
42020 - 0000	Ripette pro®	1	5

liquid handling
robotic consumables
clinical products

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- *Due to its low weight the Ripette pro is most-suitable for long dispensing series.*
- *Ergonomic design which allows an one-hand-operation of volume selection, loading and dispensing of liquids.*
- *The dosing table shows all dosing possibilities at a glance for choosing the most-suitable tip for your application.*
- *The Ripette pro is reliable and maintenance-free due to its robust design and a minimum of wear parts.*
- *By using selected premium materials the Ripette pro shows a high resistance towards a broad range of chemicals.*
- *Every device will be tested in terms of functionality and precision after production. Only after the requirements of DIN EN ISO 8655 have been fulfilled, a serial number will be issued.*

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